


**NiFTY**  
NICKEL FERRULE TECHNOLOGY

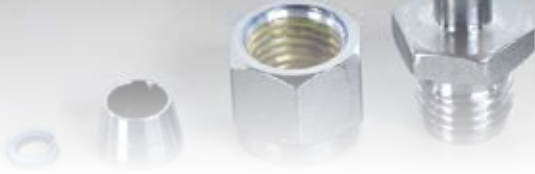

**ENTECH**  
INSTRUMENTS

## You May Notice Something New!

The ferrule, a small but critical component to leak tight sampling, has evolved. Your order is now using the latest ferrule technology from Entech.

Through years of development, Entech has created a pure Nickel ferrule that improves overall performance when compared to any other ferrule currently used for whole air sampling and analysis. The unique properties of Nickel create the best combination of inertness, low carryover, and replace-ability, so you never have to throw away a fitting, flow controller, (or worse) a 6L canister because your ferrule or ¼" Swagelok™ style nut failed you. Nickel ferrules seal and are easily replaceable just like Graphite-Vespel™ ferrules, yet just like stainless steel ferrules, they are also non-contaminating and cannot be removed by hand.




Nickel ferrules seal and are easily replaceable just like Graphite-Vespel™ ferrules, yet just like stainless steel ferrules, they are also non-contaminating and cannot be removed by hand

### ADDITIONAL PARTS

NiFTY – Nickel Ferrules (10 Pack)  
Part# 30-40400 – 1/4"

NiFTY - Ferrule Remover Part#  
30-40900 – 1/4"

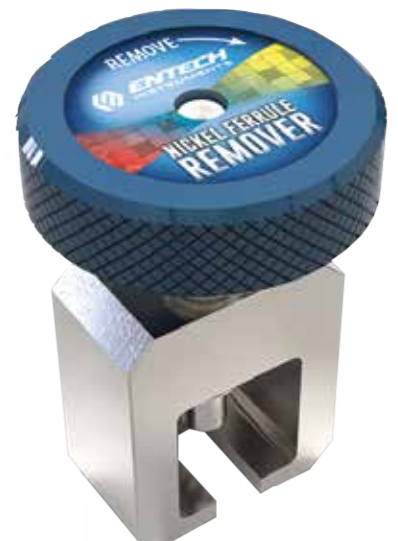

See instructions on back →

# Installing Nickel Ferrules



# Removing Nickel Ferrules

